


### Download Instructions

Right mouse click on the zip file and save it to your computer hard drive. Right mouse click again to extract the file(s), and copy the extracted file update.ini and the folder update without a subdirectory (see screenshot below) to either a CD or a USB stick.

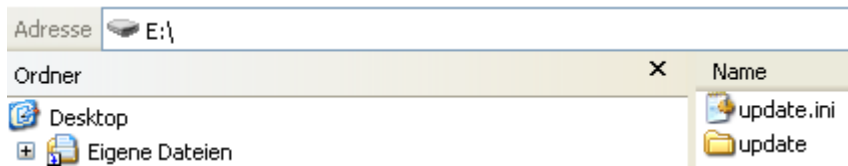

### Installation Instructions for the MagNA Pure LC 2.0 Instrument:

As new or updated purification protocols are created, they can be added to the MagNA Pure LC 2.0 Instrument software via a CD or USB stick during the Instrument start. The update media will automatically be checked for accuracy and for viruses by the Instrument software.

1. Ensure that the MagNA Pure LC 2.0 Instrument is switched off.
2. Insert the USB stick into the USB port at the front of the instrument..
3. Press the 'Power On' button. If using a CD for the update, load the CD into the CD drive as soon as the LED of the CD drive lights up.
4. The Software is automatically updated.
5. After the Main Screen is displayed, remove the update media.
6. The update is complete.
7. The Instrument is ready for use.
8. You will find the new or updated protocol in the pull-down list of protocols of the 'Ordering' sub-tab.

For details, refer to the MagNA Pure LC 2.0 Instrument Operator's Manual.